

methodbook.net

[Home](#) [Bacteria](#) [Cell culture](#) [DNA](#) [Electronics](#) [PCR](#) [Physics](#) [Probes](#) [RNA](#) [Virus](#) [Search](#)

No responsibility is assumed by methodbook.net for any injury and/or damage to persons or property as a matter of products liability, negligence or otherwise, or from any use or operation of any methods, products, instructions or ideas contained in the material herein. It is the users responsibility to ensure that all procedures are carried out according to appropriate Health and Safety requirements. Some countries have allowed a patent on the PCR process. This patent is owned by Hoffman la Roche. Make sure you have an appropriate license or are exempt before performing any PCR

Name	Institution	e-mail (optional)
Matt Lewis	Department of Pathology University of Liverpool	m.lewis@liv.ac.uk

PCR mutagenesis

Notes

This method uses a proof-reading polymerase to read all the way around a plasmid and thus incorporate the primer as the new (mutant) sequence. Only a few (say 12) PCR cycles are performed on a relatively large amount of plasmid template to minimise the chance of expanding PCR sequence errors.

Designing the primers

You need two primers, complementary to each other, containing the new (mutant) sequence flanked by 20 bases on each side. For example, suppose you have the following sequence in some gene in some plasmid;

CTA CTT CCA GAG ACA ACT GAT CTC TAC TGT TAT GAG CAA TTA AAT GAC AGC GGG

And you want to change it to;

CTA CTT CCA GAG ACA ACT GAT CTC TAC **GGT** TAT GAG CAA TTA AAT GAC AGC GGG

One primer will be;

.....5' .. CCA GAG ACA ACT GAT CTC TAC GGT TAT GAG CAA TTA AAT GAC AGC 3'

and the other primer will be the exact complement;

.....5' .. GCT GTC ATT TAA TTG CTC ATA ACC GTA GAG ATC AGT TGT CTC TGG 3'

In the PCR reaction most of these primers will be annealing to each other (and thus not extending - hence avoid normal Taq) whilst a few will be annealing to your target sequence with a small mismatch in the middle. The primers must be FPLC, HPSF or PAGE gel purified (to avoid annoying n-1 primers). Contrary to popular opinion there is no need to phosphorylate the 5' ends.

Protocol

Set up a PCR reaction like this one, use a proof-reading polymerase such as Pfu polymerase.

- 5ml 10 x Pfu polymerase buffer (inc Mg)
- 4ml 10mM dNTPs (i.e. a mixture of all four at 10mM each)
- 0.2 ml Primer 1 (primers are ~42 mers at 100pmol/ml)
- 0.2 ml Primer 2
- 1ml plasmid template (10ng)
- 37.6ml H₂O

- 2ml Pfu polymerase

The Pfu polymerase contains a 3'-5' exonuclease and it will start to chew up single-stranded DNA (the primers) from the 3' end making them shorter and less specific. Therefore assemble the reaction on ice and only add the enzyme last thing. Keep on ice until you put it into the PCR machine, already pre-heated to 94C.

PCR program

Duration	Temperature	Cycles
60 seconds	94C	1
30 seconds	94C	
30 seconds	55C	12
12 minutes	68C	

The extension temperature is 68C to minimize 'breathing' of the 5' end of primer. Then, when the polymerase has read all the way around the plasmid it is less likely to displace the primer and incorporate the 'old' sequence from the plasmid. The extension time is 2 minutes per kb of plasmid.

Cool the PCR reaction to room temperature and add 1ml of restriction enzyme DpnI.

Incubate at 37C for 1 hour.

DpnI is a 4-cutter which only cuts dam methylated DNA. The parental plasmid DNA will be cut to pieces whilst the nascent PCR DNA is left intact. All routine E.coli strains have an intact dam methylase system. Check in the back of the NEB catalogue if unsure.

Transform 5ml and 45ml of the reaction into competent E.coli.

There is no need to carry out a ligation before the transformation. The PCR product is a 'nicked' circle with the nicks in opposite strands displaced by 42bp. This is identical to a classical de-phosphorylated vector plus insert transformation (with a 42bp insert).

Miniprep some colonies, check that they are the expected size and screen 2 of them by sequencing.

I've done this twice and each time both colonies were the correct thing. I then subclone the gene out of the plasmid back into the vector in the classical way to avoid sequencing the whole vector for any PCR errors.

No responsibility is assumed by methodbook.net for any injury and/or damage to persons or property as a matter of products liability, negligence or otherwise, or from any use or operation of any methods, products, instructions or ideas contained in the material herein. It is the users responsibility to ensure that all procedures are carried out according to appropriate Health and Safety requirements.

